# HIRSCHMANN®

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## Instruction manual

Please read before use!

# pipetus®

Akkubetriebene Pipettierhilfe  
Battery operated pipette filler

Hirschmann Laborgeräte GmbH & Co. KG  
Hauptstraße 7-15 • 74246 Eberstadt Germany  
Fon +49 7134 511-0 • Fax +49 7134 511-990  
[www.hirschmannlab.com](http://www.hirschmannlab.com) • [info@hirschmannlab.com](mailto:info@hirschmannlab.com)

X2120470000

## Inhalt

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## Contents

	Page
1. Safety instructions.....	3
2. Components.....	7
3. Functional principle .....	8
4. Prohibited operations .....	10
5. Restricted operations .....	11
6. Preparation for use.....	12
7. Pipetting.....	15
7.1 Place the pipette .....	15
7.2 Fill the pipette .....	16
7.3 Empty the pipette.....	17
8. Spare parts replacement.....	19
8.1 Replacement of pipetholder complete .....	19
8.2 Replacement of filter.....	19
8.3 Replacement of safety valve.....	20
8.4 Safety valve - Important hint for assembling.....	21
9. Maintenance, cleaning and sterilisation.....	23
10. Test of tightness .....	25
11. Repair at the manufacturer .....	25
12. Trouble shooting .....	29
13. Technical data.....	31
14. Accessories – spare parts .....	33
15. Clearance certificate form.....	35

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---

### **Attention!**

Before use please read the safety instructions and observe them carefully. Before use every member of staff must be familiar with these safety guidelines and they must be kept available at all times. This instruction manual can not, however, cover every possible safety hazard. It is every user's responsibility to observe all health and safety precautions and to determine and observe restrictions before use of the unit.

- 
- Observe all general safety practices and guidelines for laboratories, such as the use of protective clothing, safety goggles and gloves when handling hazardous chemicals.
  - Never point the pipette in anyone's direction when in use!

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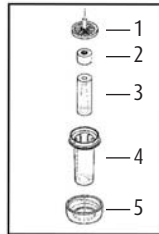
## 1. Safety instructions

- To avoid the risk of an electrical shock, never immerse either the unit or the recharger unit in water or any other liquid!
- Please ensure that the current of the power network matches with the voltage on the charger unit.
- In the event of any damage of the cable or the recharger unit, the recharger unit should not be used.
- Never touch the plug of the recharger unit with wet hands!
- Observe the guidelines and procedures of the chemical manufacturers.
- The recharger unit may only be used in rooms. Protect from moisture!
- Non-rechargeable batteries should not be connected to the recharger unit!

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## 1. Safety instructions

- A safety valve (2) in the pipetholder case (4) is located on top of the silicon adaptor (3) which prevents the liquid from getting into the handle, if the pipette is accidentally overfilled. If the safety valve has been activated, it must be removed by unscrewing the pipetholder and then cleaned and dried. Afterwards it must be inserted in the prescribed direction (see point 8). Wear safety gloves and goggles! **(Danger of infection)**

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---

### **Attention!**

If the unit is opened, the warranty is invalidated and any resultant claims void.

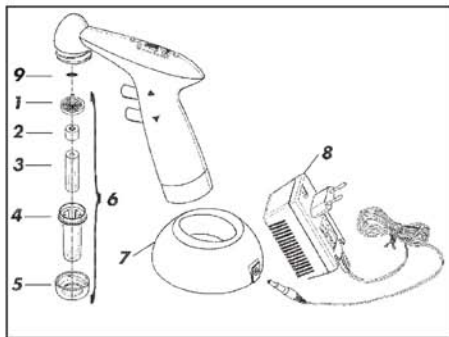
---

- The filter supplied contains a hydrophobic 0,2 µm membrane for air filtration in either direction.
  - Only have the unit professionally serviced.
  - Never use undue force!
- 

### **Attention!**

Only ever use the original recharger unit, accessories and replacement parts to avoid damaging the unit.

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
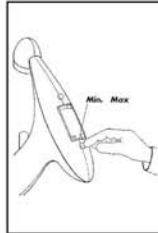
### 3. Functional principle

- The unit is a rechargeable pipetting controller for use with all standard pipettes from 0,1 ml to 200 ml.
- The pump is activated by pressing the In-( $\triangle$ ) or respectively the Out-( $\nabla$ ) button. The speed of filling (In) or dispensing (Out) is regulated by the intensity of pressure on the buttons.
- The pipetting speed is set by switching the 5 speed-stages and therefore an individual adaption to the pipetted volume is possible.

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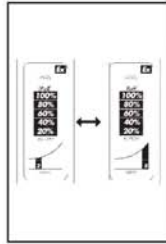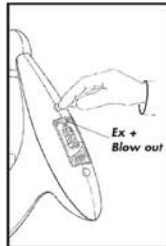
### 3. Functional principle

- The selected speed stage is indicated at the bottom of the LC display.

- By switching from “Ex” to “Blow out” or vice versa, the unit can accommodate both to deliver and blow-out pipettes. In the “Ex” position the pump remains switched off while atmospheric pressure empties the pipette. In this way, the working time of the batteries is doubled.

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---

##### **The unit can not be used with:**

- liquids of which vapours attack the materials silicone, PTFE, Novodur, PVDF, NBR, PCPA, PP and PA6.
  - easily inflammable liquids (Attend to the flame point).
  - usage in explosive areas.
-

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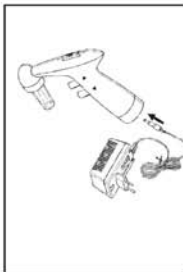
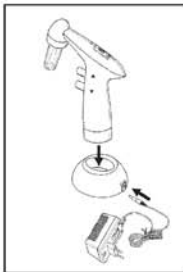
## 5. Restricted operations

- Working and charging temperatures between +10 °C and +40 °C.
- The unit should only be used and optional batteries loaded in areas where it can be assured that no hazardous potentially-explosive atmospheres occur.
- Load only in Ex-secure areas.
- Do not keep unsupervised in an explosive area.

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

## 6. Preparation for use

Before putting the unit into operation, the batteries must be charged:

- Either through inductive charging technology in the charging stand (time for charging approx.  $\leq 8$  h). Insert the plug of the charger unit into the socket of the charging stand.
- Or through direct charging of the unit (time for charging approx.  $\leq 8$  h). Insert the plug of the charger unit directly into the socket of the unit.
- Insert the plug of the charger unit into a table or wall socket.

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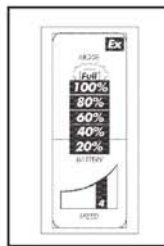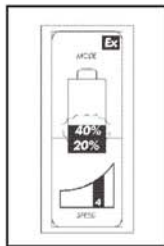## 6. Preparation for use

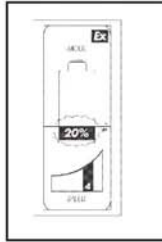
- One after the other segment of the battery capacity indication starts flashing during the charging procedure according to the respective charging status. Provided e. g. that the batteries are almost empty, the segment 20 % will be flashing.
- If the batteries are charged completely the display "Full" will be flashing.

### Guide:

The warming up of the unit during the charging operation is harmless.

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#### 6. Preparation for use

- During pipetting one after the other segment of the display will disappear.
- When falling short of 10 % of the battery capacity, an acoustic signal will be heard after the pipetting operation.
- This indicates, that the unit must be charged again as soon as possible.

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---

### **Attention!**

The unit may only be charged with the original recharger unit. The network voltage must match with the voltage indicated on the recharger unit (maximum deviation  $\pm 10\%$ ).

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## **7. Pipetting**

### **7.1 Place the pipette**

- Place the pipette into the holder to ensure a secure and airtight fit!

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---

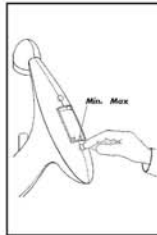
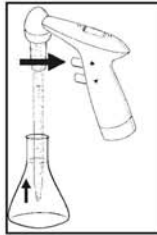
### **Attention!**

Danger of glass breakage, especially with thin pipettes.

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## 7.2 Fill the pipette

- To fill the pipette, press the In-button, the intensity of finger pressure on the button regulates the working speed.
- Select one of the 5 speed stages according to the volume of the pipette.

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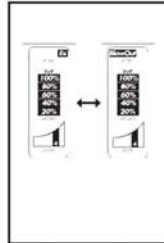
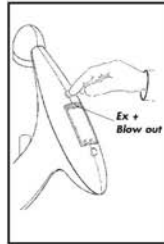
### **Attention!**

No liquid should get into the unit as this can adversely affect the operating functions.

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
### 7.3 Empty the pipette

- If Blow out pipettes are used, the key has to be set so that “Blow out” appears in the display.
- If To Deliver pipettes are used, the key has to be set so that “Ex” appears in the display.
- The chosen pipetting mode is indicated in the upper part of the display.

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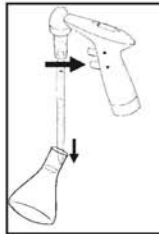### 7.3 Empty the pipette

- => To empty the pipette, press the  
▼ button, the intensity of finger pressure on the button regulates the working speed.

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---

#### Guide:

- Pipetting mode Blow out => pipette is blown out by the pump.
  - Pipetting mode Ex => pipette is emptied by atmospheric pressure (without the pump).
-

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## 8. Spare parts replacement

### 8.1 Replacement of pipetholder complete

- Screw off the nut.
- Replace the pipetholder by a new one.

### 8.2 Replacement of filter

- Screw off the nut.  
Remove the old filter and place the new one in the pipetholder case so that the pivot shows upwards. Screw the pipetholder onto the unit, by turning the nut clockwise.

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### 8.3 Replacement of safety valve

- Screw off the nut (5) counter-clockwise. Take out the filter (1) and place it aside. After that, remove the old safety valve (2) and place the new one in the pipetholder case (4), **so that the black sealing ring shows upwards**. Place filter (1) on the safety valve with the pivot facing up. Screw the pipetholder onto the unit, by turning the nut clockwise.

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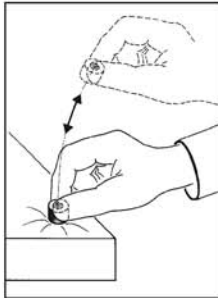
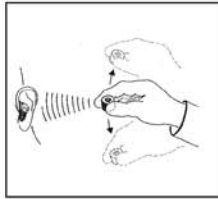
## 8.4 Safety valve

### IMPORTANT HINT FOR ASSEMBLING

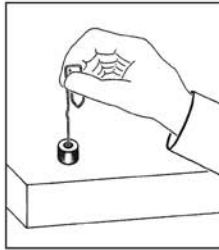
Black O-Ring of safety valve must be visible upside above silicone adapter!

Function check of the safety valve. The membrane of the valve must create a “rattling” sound, while shaking the valve.

A sticking valve might become flexible again, by gently knocking it e. g. on the table edge. However in the event that the valve is glued, it must be soaked in a cleaning solution, rinsed with water and dried afterwards. Do not open the valve.

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
#### 8.4 Safety valve

In case the function of the valve is not restored yet, try to make it work again, by gently pushing the membrane with a thin and blunt object.

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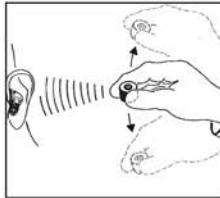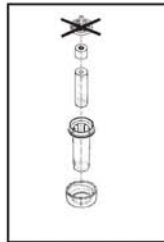
## 9. Maintenance, Cleaning and Sterilisation

- Under normal conditions of use, the unit is maintenance-free. In the event, nevertheless, that they do become dirty or the suction becomes impaired, the safety valve and membrane filter should be replaced and the unit recharged.
  
- The pipetholder, except the membrane filter, can be sterilised at 121 °C.

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- The safety valve (2) must be cleaned and dried after autoclaving. While reassembling the unit make sure that the **side of the safety valve (2) with the black sealing ring shows towards to the filter (1)**.

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---

### Attention!

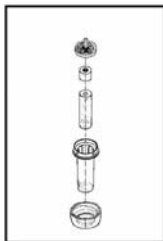
When cleaning and exchanging the safety valve and membrane filter, protective gloves and goggles must be worn (Danger of infection)!

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## 10. Test of tightness

Once the required volume has been aspirated into the pipette, the liquid has to be kept in it unless the buttons are pressed. If the pipette however is dripping, disassemble the pipette holder and reassemble it again carefully.

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## 11. Repair at the manufacturer

Please observe the following points should it prove necessary to return the unit for repair (despite having handled it with all due care):

- Before sending the unit back, the user must ensure that it is carefully cleaned and decontaminated. The healthwise harmlessness must be confirmed, send the confirmation together with the unit to the manufacturer or the dealer, (a copy of the clearance certificate can be found on page 35).

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#### 11. Repair at the manufacturer

- Units will only be repaired, when they have been cleaned and decontaminated by the user.
- For safety reasons dirty units will be returned to the sender without being repaired.
- Please describe the malfunction exactly.
- Please quote the pipetted liquid.
- In the event of the unit being tampered with or being used in any improper way, (e.g. the application of undue force) any claims under the warranty are void.

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## 12. Trouble shooting

Problem	Possible reason	Solutions
Liquid column in pipette not maintained (liquid drips or runs out of pipette)	<ul style="list-style-type: none"><li>- Safety valve is incorrectly inserted</li><li>- Pipette holder leaks</li><li>- Seal ring in housing (on transition from filter to housing) is defective, incorrectly inserted or missing</li></ul>	<ul style="list-style-type: none"><li>- Insert the safety valve so that the side with the black ring (O-ring) faces in the direction of the filter (see 8 and 9).</li><li>- Check the sequence for pipette holder assembly. Check the function of individual parts where appropriate and replace if necessary.</li><li>- Stand the unit on its head and check the seal ring in the housing above the pipette holder to ensure it is fitted correctly, or insert if necessary.</li></ul>
The pump will not blow out the pipette	<ul style="list-style-type: none"><li>- Incorrect pipetting mode</li></ul>	<ul style="list-style-type: none"><li>- Set the Ex/Blow out key to the Blow out mode. The pipette will be blown out!</li></ul>
Note: Pipetting mode on Blow out	=> the pump will blow out the pipette	
Pipetting mode on Ex	=> the pipette runs out via atmospheric pressure (without pump)	

## 12. Trouble shooting

Problem	Possible reason	Solutions
The pump does not function neither filling nor blowing out	<ul style="list-style-type: none"><li>- Batteries flat</li><li>- Safety valve incorrectly positioned</li></ul>	<ul style="list-style-type: none"><li>- Recharge the batteries with the re-charger unit.</li><li>- Place the safety valve so that the black sealing ring shows towards the filter.</li></ul>
Overfilled pipette	<ul style="list-style-type: none"><li>- Safety valve is activated immediately and prevents liquid from getting into the unit</li></ul>	<ul style="list-style-type: none"><li>- Remove safety valve, clean and re-assemble. If required, replace the safety valve. During assembly take note of point 8.3.</li></ul>
Suction too slow or too weak	<ul style="list-style-type: none"><li>- Selected speed stage is too low</li><li>- No tight, secure connection of the pipetholder</li><li>- Filter is full</li><li>- Silicone adaptor is damaged</li></ul>	<ul style="list-style-type: none"><li>- Improve the speed selection to a greater stage e.g. 3-&gt;5.</li><li>- Screw the pipetholder more tightly onto the handle.</li><li>- Exchange the filter.</li><li>- Exchange the silicone adaptor.</li></ul>

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#### Batteries:

- 2 NiMH batteries
- Operation time for permanent use:
  - Position Ex 50 h
  - Position Blow out 25 h
- time for charging:
  - inductive charging technology  $\leq 8$  h
  - direct charging  $\leq 8$  h

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### 13. Technical data

CE:

This mark in the confirmation that the unit conforms to the established requirements of the EU guidelines and has been tested according to the criteria.

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## 14. Accessories – spare parts

Silicone adapter	990 32 01
Safety valve	990 32 02
Pipetholder complete	990 38 00
Pipetholder case	990 38 03
Nut	990 38 02
Round filter set (consisting of 5 replacement filters)	990 37 04
Recharger unit complete with inductive Charging stand, 100-240 V/50 Hz	990 75 03
Stand	990 57 00
O-ring	990 32 05

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## 15. Clearance certificate form

for unit ..... (Serial number) Medium used: .....

	YES	NO	Please define in case of Yes
Blood, bodily fluids, patholog. samples:	<input type="checkbox"/>	<input type="checkbox"/>	_____
Other hazardous biologic substances:	<input type="checkbox"/>	<input type="checkbox"/>	_____
Biologically-degradable materials which can pose a risk:	<input type="checkbox"/>	<input type="checkbox"/>	_____
Chemicals containing substances harmful to health:	<input type="checkbox"/>	<input type="checkbox"/>	_____
Radioactive materials:	<input type="checkbox"/>	<input type="checkbox"/>	_____
Other hazardous and toxic substances:	<input type="checkbox"/>	<input type="checkbox"/>	_____

### We hereby declare that

- the unit has been carefully cleaned and decontaminated prior to shipment
- no risks resulting from contamination exist
- we are aware that liability claims can be made against us in the event of damage being caused by contaminated units.

### Address of sender:

Name: \_\_\_\_\_

Date: \_\_\_\_\_

Signature: \_\_\_\_\_